RANGER[®] 7000
Ethernet Interface
Instruction Manual

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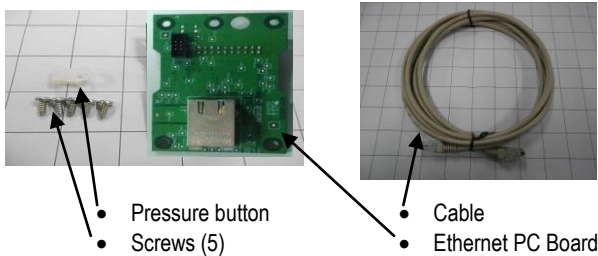
INTRODUCTION

This Ethernet interface kit is for use with the Ranger[®] 7000 scales.


KIT CONTENTS


INTERFACE INSTALLATION

Only one interface module can be installed in the terminal option slot at a given time. Install the Ethernet Interface on the underside of the balance as shown.

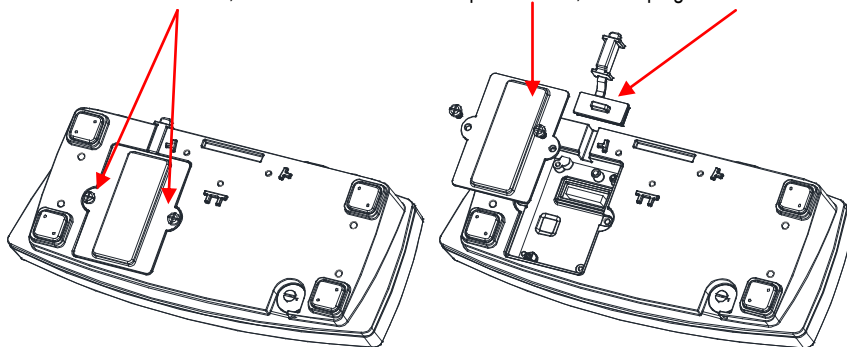1. Release the two screws, remove the red interface options cover, and unplug the black rubber insert.

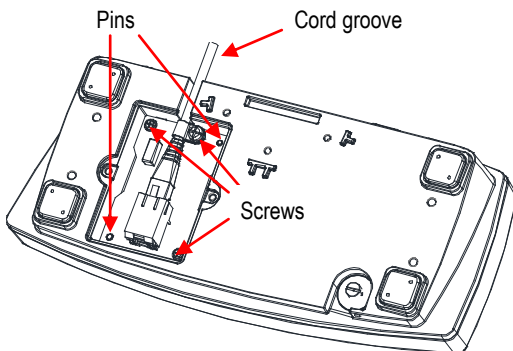2. Install the Interface PCB. Push it through the pins and secure it with the 3 screws.

3. Connect the Ethernet Cable.
4. Place the cord in the groove (see step 2) and put the red interface option cover back on.

SETUP

Upon installation, the scale will recognize the Ethernet Interface Kit and add relevant items to the menu. Configure the scale to the desired RS232 and printing parameters (refer to the scale User Manual for assistance in using the menus).

Configuration	
Host Name	Host Name
MAC Address	00:E0:7C:00:00:00
Port	9761
Version	1.00
DHCP	On >
IP Address	169.254.1.1

Print Setup

	Reset	>
	Stable Weight Only	Off >
	Auto Print	>
	Print Template	Simple >
	Edit Template	Simple >
	Line Feed	1 Line >

Configuration	
IP Address	169.254.1.1
Gateway	169.254.1.1
Subnet Mask	255.255.255.0
Primary DNS	169.254.0.1
Secondary DNS	169.254.0.2

Ethernet Connection

If DHCP is set to ON, the IP Address will be automatically assigned by the IP share or router. If the DHCP is set to OFF, the user can setup the IP Address, Subnet mask, Gateway, Preferred DNS and Alternate DNS. The Interface has a preset IP Address. The Port Number is always **9761** and cannot be changed.

Note: After the setup of IP Address, Gateway, Preferred DNS and Alternate DNS, please reboot the scale to enable the setting.

OPERATION

Output

With the interface installed the terminal will operate in one of the following two ways:

- Auto Print = off, Stable = on or off
Pressing PRINT will send the display data.
If Stable is set to On, the balance will wait for a stable reading before sending the stable data.
- Auto Print = not off, Stable = on or off
The balance will automatically send data based on the settings in the Menu.
If Stable is set to On, only stable values will be sent.

Data sent from the Interface is in standard ASCII format terminated with a [CRLF] (carriage return-line feed). For the output format please refer to the Ranger 7000 instruction manual.

Input

The scale will respond to various commands sent via the interface adapter. Terminate the following commands with [CRLF].

Commands

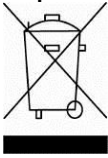
For the full list of commands, please refer to the Ranger 7000 instruction manual.

ACCESSORIES

For a complete listing of OHAUS printers and other accessories, contact OHAUS Corporation or visit www.OHAUS.com

COMPLIANCE

Disposal

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to www.ohaus.com/weee.

Thank you for your contribution to environmental protection.

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