

FLEXIBLE-ARM™ LIGHT INSTRUCTIONS

SPECIFICATIONS:

Electrical: 150 VAC, 60Hz, UL Listed
 Light: 35 watt Halogen bulb
 Plug: Polarized
 Arm length: **Cat. No. 3095** - 21"
 Cat No. 3096 - 31"
 Weight: 8 lbs.

FEATURES

- Base - 7" diameter, 7lb.
- 35 watt bright Halogen white light
- Clamping unit for vertical or horizontal Positioning
- Switch in cord
- Flexible-Arm™ articulates at 3 points
- Fingertip bar for raising, lowering and pivoting with ease
- Spring counterbalance for stability
- Vented shade with rolled edges
- Safety-glass with UV filter

BASE ASSEMBLY (see Figure 1)

1. Remove "Screw I" from base "Stem II".
2. Insert the lamp frame into base "Stem II", align screw holes.
3. Screw in "Screw I" and finger tighten.

CLAMPING UNIT ASSEMBLY

1. Insert the lamp frame into hole in top of clamp assembly.
2. Loosen clamp with lever on bottom of assembly.
3. Position clamping mechanism and tighten firmly using lever on bottom of assembly.

BULB REPLACEMENT (see Figure 2)

1. Turn off lamp and allow shade and accessories to cool.
2. Turn the lampshade upward.
3. Remove CENTER screw (using a Phillips screwdriver) and clip from protective glass cover.
4. Pull out failed bulb and replace with halogen bulb. (Control Company Cat. No. 3098)
5. Replace protective glass and clip and screw in place.

NOTE: Do not remove screws holding fingertip bar in place.

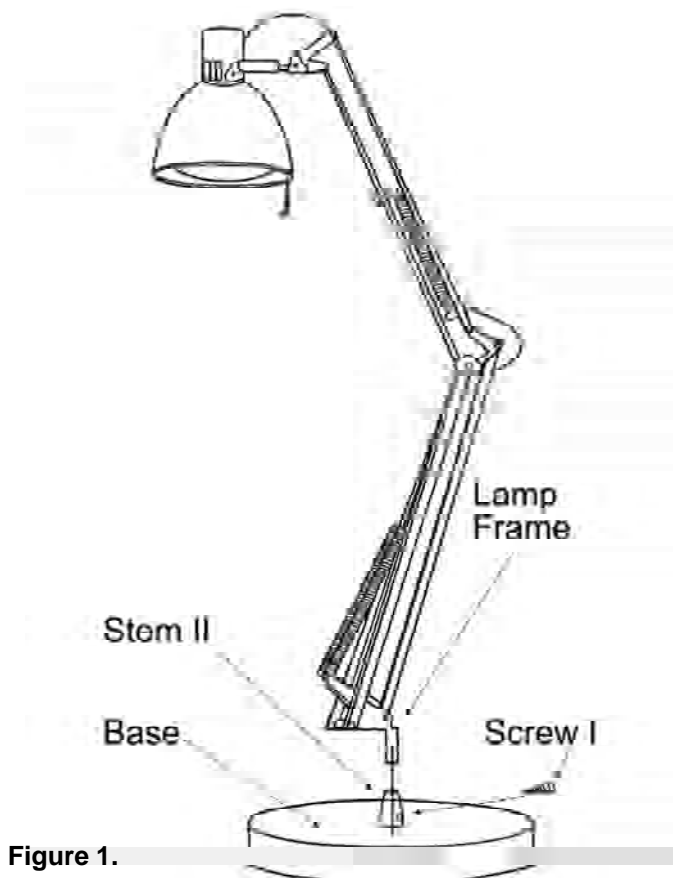


Figure 1.

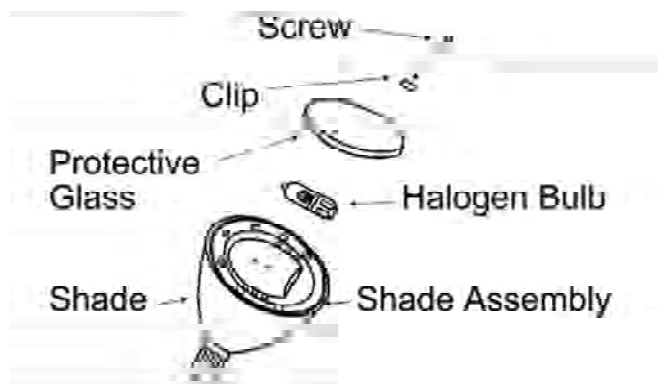


Figure 2.

CONTROL COMPANY

Control Company is
 ISO 9001 Quality-Certified by DNV and
 ISO 17025 accredited as a
 Calibration Laboratory by A2LA.

Cat. No. 3095/3096
Flexible-Arm™ Light Instructions
92-3095-10
 Rev. 1 071504

Traceable® is a registered trademark of Control Company



© 2004 Control Company

1.888.610.7664



www.calcert.com

sales@calcert.com