

Traceable®

4-Channel Timer Quick-Start Guide



WARRANTY, SERVICE, OR RECALIBRATION

Traceable®
an antylia scientific company


Traceable® Products are ISO 9001:2018 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Item No. 56000-16

Model No. 5667

©2022 1065T9_M_92-5667-00 Rev. 0 09022022


SETTING A TIMER

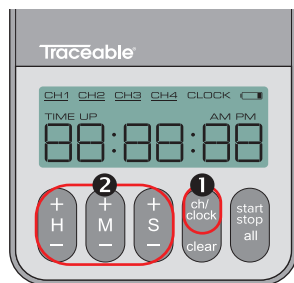

1. Press CH/CLOCK button until the desired channel is solid (not flashing) on the screen.
2. Use the H/M/S buttons to configure timer.


Note: To clear and start over, press CLEAR button.

STARTING A TIMER

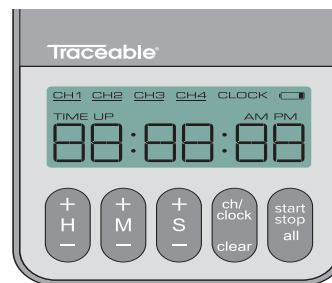


- When starting/stopping an individual timer, navigate to that timing channel using the CH/CLOCK button and then press START/STOP ALL.
- To start/stop all timers, press the START/STOP ALL button from the clock screen.
- Timers which are counting down will lash quickly.
- Timers which are counting up will lash slowly.
- Timers which have expired will lash rapidly.

Traceable® 4-Channel Timer

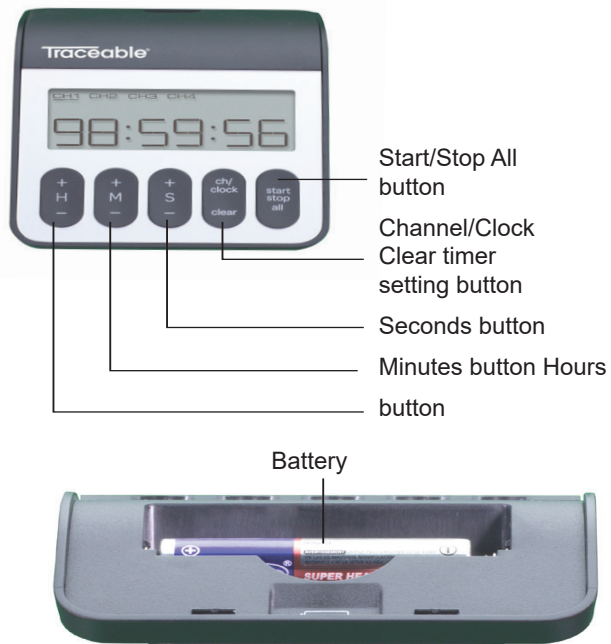

INITIAL SETUP

- 1 Remove battery cover and insert batteries

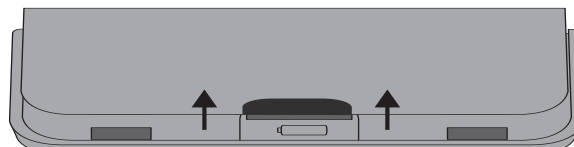

- 2 Replace battery cover
- 3 Press CH/CLOCK button until clock appears on top row of display.



Hold CH/CLOCK until there is an audible beep and the time flashes



- 5** To adjust 12/24 hour clock, press and hold START/STOP ALL button.



- 6** Use H/M/S buttons to adjust the time



- 7** Pressing Seconds button will round to the nearest minute.

- 8** Press the CH/CLOCK button to exit setup.